



Techne qPCR lyophilised 2X qPCR MasterMix

Instructions for use of Techne Prime Pro lyophilised MasterMix

Contents

Introduction	3
Kit Contents	4
Kit Storage	4
Suitable Sample Material	4
Licensing Agreement and Limitations of Use	4
Techne Satisfaction Guarantee	4
Quality Control	4
Bench-side Protocol	5
Reaction setup	6
Amplification Protocol	7

Introduction

Techne Prime Pro 2X lyophilised MasterMix is optimised for use in qPCR. The product is stable at ambient temperatures for at least 18 months and can be conveniently shipped and stored at room temperature. The MasterMix contains a thermo-stable TAQ Polymerase as well as buffer, MgCl₂ and stabilisers at concentrations optimised for the enzyme. The MasterMix has been freeze-dried to produce a room temperature stable preparation.

The kit includes the lyophilised MasterMix, re-suspension buffer and a tube of ROX dye which can be added as required when the MasterMix is to be used on hardware platforms that use ROX as a passive reference dye.

The re-suspended solution requires only the addition of your cDNA, primers and probe to be PCR ready.

The performance of Prime Pro lyophilised MasterMix is as good as or better than leading brands.

The perfect partner for Prime Pro kits

Techne Prime Pro 2X lyophilised MasterMix is designed for use with our range of Prime Pro qPCR pathogen detection kits. Together they represent the ultimate solution for convenient logistics as well as high quality qPCR performance.

Kit Contents

- 3 x Lyophilised MasterMix (50 reactions per glass ampule)
- 1 x Lyophilised ROX (**BROWN**)
- 4 x Re-suspension buffer (**BLUE**)

Kit Storage

The Prime Pro Lyophilised MasterMix should be stored at ambient temperature on arrival. The kit is stable for at least 18 months at ambient temperature. Once resuspended in the provided buffer the kit should be stored at -20°C. Repeated freeze/thawing will not compromise the performance of the product. Under these conditions reagents are stable for six months from date of resuspension.

Suitable Sample Material

All kinds of sample material suited for PCR amplification can be used. Please ensure the samples are suitable in terms of purity, concentration and DNA integrity. Always run at least one negative control with the samples. To prepare a negative-control, replace the test sample with RNase/DNase free water.

Licensing Agreement and Limitations of Use

PCR is covered by several patents owned by Hoffman-Roche Inc. and Hoffman-LaRoche, Ltd. Purchase of Techne kits does not include or provide licence with respect to any patents owned by Hoffman-La Roche or others.

Techne Satisfaction Guarantee

Techne takes pride in the quality of all our products. Should this product fail to perform satisfactorily when used according to the protocols in this manual, Techne will replace the item free of charge.

Quality Control

As part of our routine quality assurance programme, all Techne products are monitored to ensure the highest levels of performance and reliability.

Bench-side Protocol

1. For each glass ampule; Re-suspend lyophilised MasterMix in 525µl of re-suspension buffer

Do not replace the re-suspension buffer with water or any other buffer.

The MasterMix is then ready to use as a 2X qPCR MasterMix

2. Add ROX if required

ROX is required for platforms that use ROX as a passive reference guide. Use table 1 below to see if ROX addition is required for your hardware platform. If ROX is required then follow the instructions below.

- Re-suspend the Lyophilised ROX (**BROWN**) in the correct volume of re-suspension buffer (**BLUE**) according to table 1 below.
- Add re-suspended ROX to each ampule at the correct level.

Table 1. ROX addition

Real time PCR platform	ROX re-suspension volume	ROX addition per ampule
Applied Biosystems 7700, 7000, and 7900, 7300 StepOne, StepOnePLUS and ViiA7 platforms, Roche capillary Lightcyclers.	100µl	20µl
All Stratagene platforms	200µl	15µl
Applied Biosystems 7500 platform	700µl	10µl
All Other machines	NOT REQUIRED	NOT REQUIRED

Reaction setup

- **When using Techne Prime Pro pathogen detection kits.**

For each 20µl real-time PCR reaction add the following to each reaction tube

Components	1 Reaction	
Prime Pro Lyophilised 2X qPCR	10	µl
Primer/Probe mix	1	µl
Template (25ng)	5	µl
RNAse/DNAse free water	4	µl
Final volume	20	µl

- **Suggested use with user supplied primers and probe.**

For each 20µl real-time PCR reaction add the following to each reaction tube

Components	1 Reaction	
Prime Pro Lyophilised 2X qPCR	10	µl
Primers (6pmols Forward and Reverse)	x	µl
Probe (3pmols)	x	µl
Template (25ng)	x	µl
RNAse/DNAse free water (up to Final volume)	x	µl
Final volume	20	µl

*6pmols of primer gives a working concentration of 300nM in a 20µl reaction

Amplification Protocol

- For use with Prime Pro pathogen detection kits

	Step	Time	Temp
	Enzyme Activation - Hot Start	2 min	95°C
Cycling x50	Denaturation	10s	95°C
	DATA COLLECTION*	60s	60°C

*Fluorogenic data should be collected during this step according to the detection kit.